



The Experts in
Medical Refrigeration

USER GUIDES

How to change between
air or load temperatures
using the controller



Ebrington
MEDICAL

1 UNLOCK THE CONTROLLER



1 Press and hold the navigation button for 3 seconds



2 Hold until all dashes appear and the screen below appears

2 CHANGING BETWEEN AIR OR LOAD TEMPERATURES



1 You can now use the $\frac{\text{AIR}}{\text{LOAD}}$ button to switch between air and load temperatures



AIR TEMP



LOAD TEMP

2 The controller will display A or L depending on the selection